


## TOCSY 2D spectroscopy

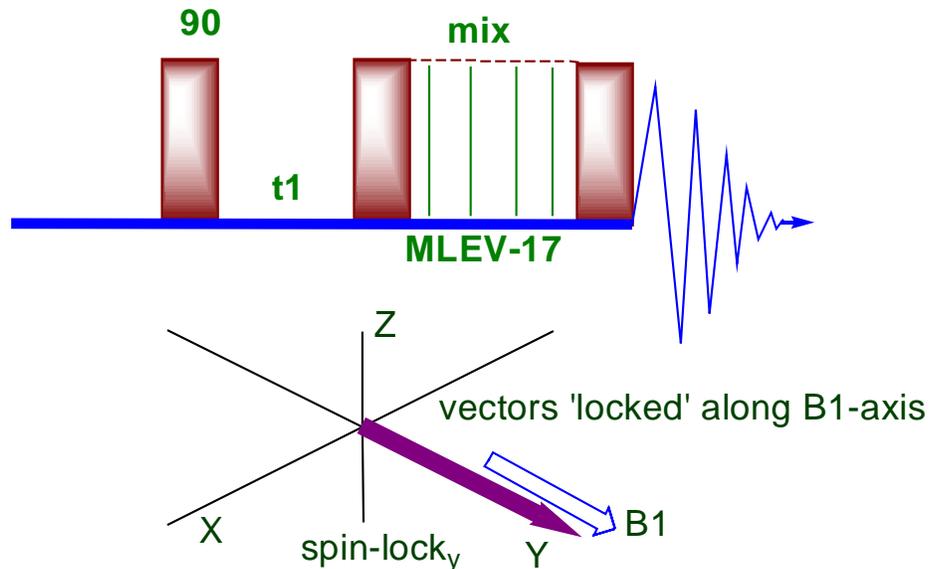

### Introduction

**TOCSY** is the abbreviation for **T**Otal **C**orrelation **S**pectroscop**Y**. It is also called **HOHAHA** (**H**Omonuclear **H**Artmann **H**Ahn). During this pulse sequence, after the evolution period  $t_1$ , the magnetization is spin-locked (for example by a series of  $180^\circ$  pulses) for a period of time on the order of 20-100 ms. During this mixing time the magnetization exchange through scalar coupling. During this spin-lock period, the magnetization behaves as a strongly coupled spin system and evolves under the influence of a "collective spin-mode". In that collective mode, coherence transfer is possible between all coupled nuclei in a spin system, (even if the are not directly coupled). One of the most popular mixing schemes in TOCSY is the MLEV-17.

For small mixing period, (e.g. 20-30 msec) Relay1-COSY type of data can be observed. As the mixing period gets longer, correlation with more distant protons can be observed (e.g. mix=80-100 msec can correlate H1 to H6 in carbohydrate). The extent of correlation depends mainly on the length of the mixing period.

This sequence is much more useful than the Relay-COSY as it very efficiently transfer magnetization before relaxation takes place.

### TOCSY experiment

1. Acquire a 1D  $^1\text{H}$  NMR first, calibrate **pw90** at **tpwr=57**.
2. Set cursors  $\sim 0.5$  ppm beyond the peaks on both sides.
3. **movesw** ↵
4. re-acquire a 1D NMR.

5. mf(1,3) jexp3 ↵
6. TOCSY ↵
7. dg ↵

You should see that **tpwr=57** with **pw=23.25**. the pw was calibrated previously at tpwr=57. the spin lock power, *slpwr*, is usually lower (46-49 dB). This experiment, it is set to 48 which is 9dB smaller than *tpwr*. So *slpw* should be increased in this way.  $23.25 * 2 * 1.414 = 65.188$ . The math is like that, if you lower 1dB, you need to increase


pulse width by  $\sqrt[6]{2} = 1.12$  times. So you have 9 dB decrease, then you need to

increase pulse width by  $(\sqrt[6]{2})^9 = 2 * \sqrt{2} = 2 * 1.414$

8. dps ↵


ACQUISITION		TRANSMITTER		FLAGS		PRESATURATION		
seqfil	TOCSY	tn	H1	hs		n	satmode	n
sw	3870.0	sfrq	399.729	sspul		y	satfrq	0
at	0.265	tof	-403.2	PFGflg		y	satdly	0
np	2048	tpwr	57	hsglvl		4854	satpwr	0
ss	64	pw	23.250	hsgt	0.002000		PROCESSING	
d1	0.600	DECOUPLER		SAMPLE		gf	0.061	
nt	2	dn	H1	date	Mar 3 2004	gfs	not used	
ct	2	dm	nnn	solvent	CDC13	fn	4096	
2D ACQUISITION		SPIN LOCK		sample	undefined	2D PROCESSING		
sw1	3870.0	slpwr	48	SPECIAL		gf1	0.048	
ni	200	slpw	65.188	temp	not used	gfs1	not used	
phase	arrayed	mix	0.080	spin	not used	proc1	lp	
		trim	0.0020	gain		fn1	2048	

9. **np=2048**, number of points in t2, it must be even number. A number to the power of 2 is required. This will also give a value to **at**. **Bigger np gives better resolution in F2 dimension.**

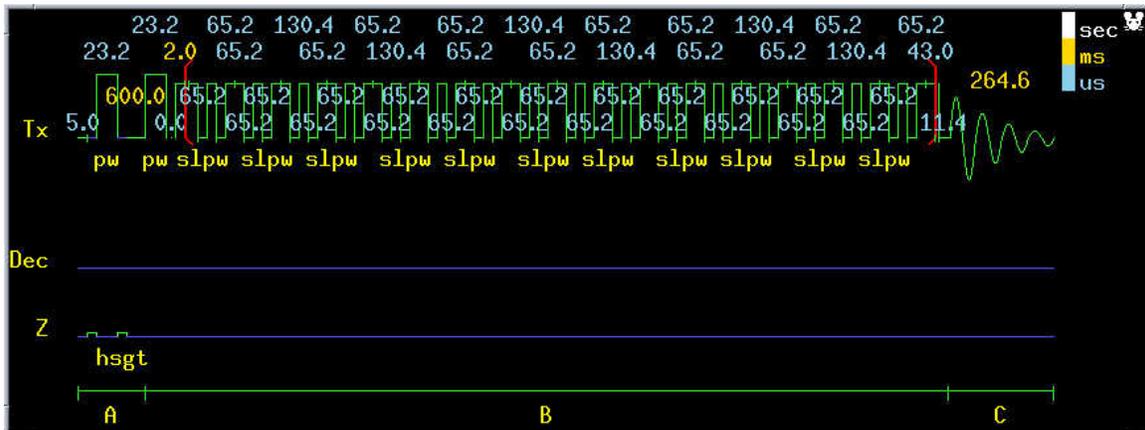


10. **phase =1,2** which is *arrayed* as is seen in dg window. Type **da** to see the details.

10.  $at=np/sw=2048/3870$ (which has been already calculated by the computer)
11. **slpw** is the spin lock pulse width which must be a 90 degree pulse width.
12. **slpwr** is the power of the spin lock, the max is 49. It must be correlated with **slpw**.
13. **mix** is the mixing time; the larger the mixing time the more transfers will occur(*the more protons will be correlated*). It is normally 0.080-0.100 s, which is suitable for 5-8 transfers. Mix is
14. **d1** can be longer than gCOSY. {**T1=200ms** for strychnine, so **d1=0.6** is OK. Normally you should set  $d1+at \geq 3T1$ . Here  $0.6+0.265 = 0.865$  which is more than  $3*T1(=0.6)$ }.
15. **ni=200** or shorter if you do not have enough time. To have a reasonable value of ni, please remember this:  $ni = sw1/12 \sim sw1/24$ . In this experiment, it is between 161~322. So 200 is a suitable value. { $sw1/(2*ni)$  is the resolution in F1, an idea value is 6Hz/pt if time allows. But it should be smaller than 12Hz/pt}.
16. turn **spin** off, set **gain** to the highest value which does not cause overflow. **go** ↓ if ADC overflow, **aa** ↓ lower gain, **go** ↓. { you can type **set2dgain** or type **ni=1 nt=1 ss=1 phase=1 gain='n' ga** then after the experiment finishes, type **gain='y'**}
17. **svf** ↓ TOCSY\_strychnine(or whatever you like) ↓
18. **setLP1 sqcosine fn=np fn1=fn wft2da** ↓

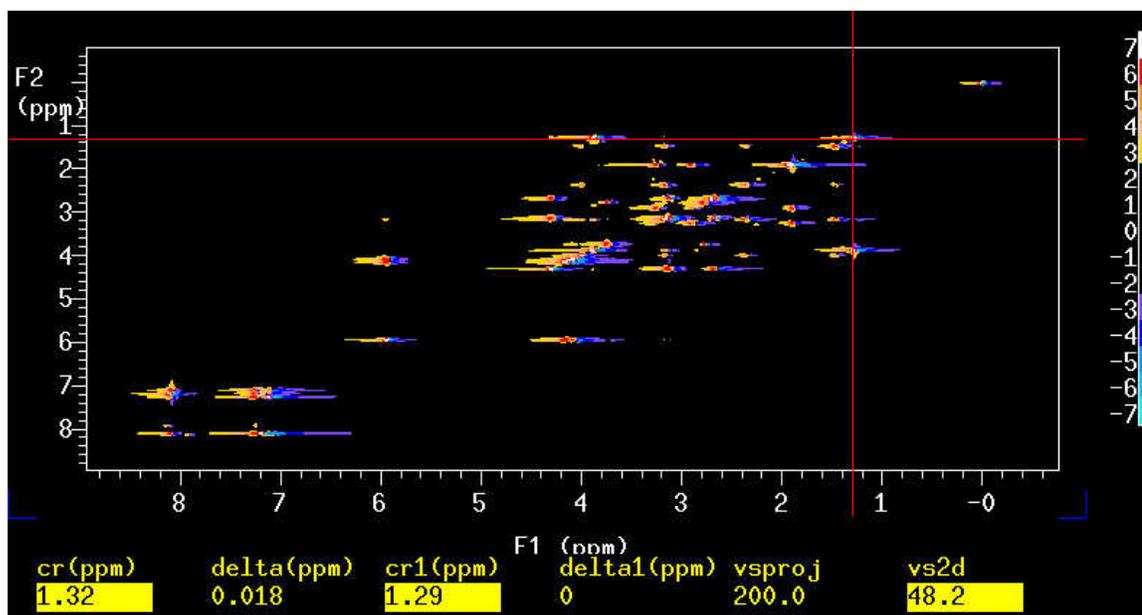


*You see the above spectrum. Apparently you need to phase the 2D spectrum.*


19. **pmode?** **pmode='full'** ↓ if it was 'partial'.
20. **trace='f1'** if it was 'f2'.
21. point the mouse at the peak of the molecule( do not point at solvent peak)
22. **ds** ↓

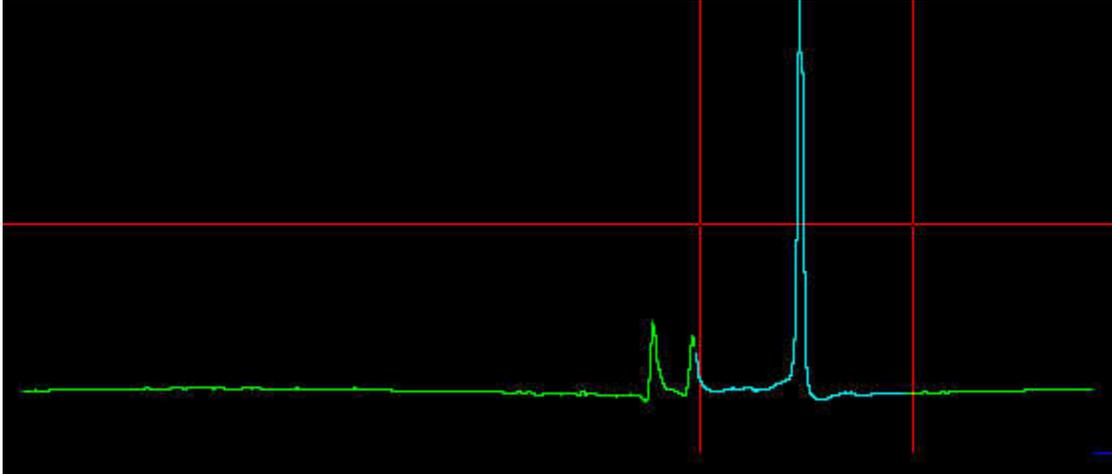

23. click phase button and phase the spectrum as you do on a normal 1D. This time you are adjusting rp1 value, the right phase on f1.
24. **dconi** ↵ point the mouse at the left peak **ds** ↵

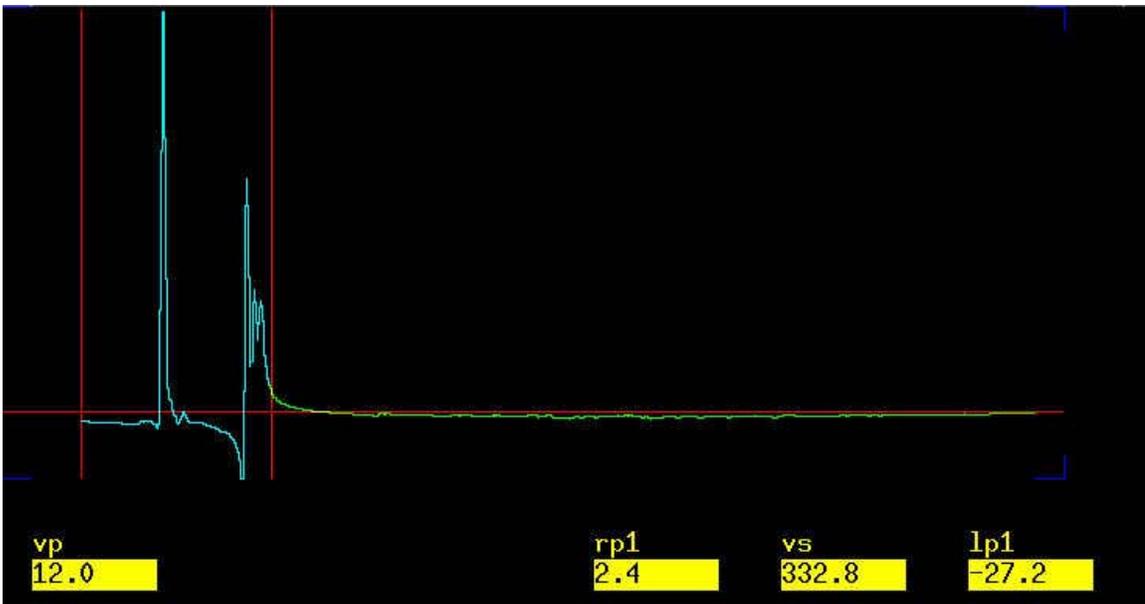

25. adjust the phase of the right side peak as shown above.
26. **dconi** ↵
27. **trace='f2'**, do the same thing as you did on f2 dimension.
28. **dconi** ↵ adjust vs **.vs+20%; vs-20%**.

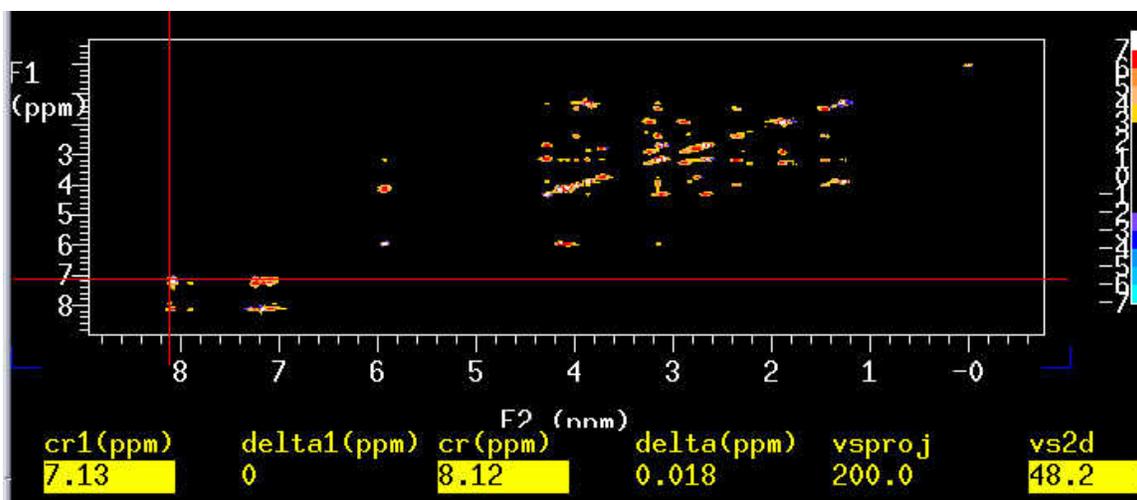

You should see the 2D like this.

29. *Re-referencing the spectrum.* Point the mouse at the CDC13 diagonal peak and type `rl(7.27p) rl1(7.27p)`.

30. `pcon(10,1.2) page`↵

31. Print with 1D on side with 1D in experiment 1: `plcosy(10,1.2,1)`.